

SOP Title: KAPA Library Quantification (Illumina Platforms)

1.0 Purpose and scope

This standard operating procedure (SOP) is for the quantification of next generation sequencing (NGS) libraries for Illumina platforms on the Applied Biosystems (ABI) ViiA7 instrument using QuantStudio Real-Time PCR Software v1.2.

Related documents:

- Manufacturer supplied protocol for library quantification: KAPA Library Quantification Kit Illumina Platforms (KR0405-v9.17). [Link to document.](#)
- Manufacturer supplied reference material: KAPA Library Quantification Technical Guide-KAPA library Quantification Kits for Illumina Platforms V1.14. [Link to document.](#)
- Manufacturer supplied: QuantStudio Real-Time PCR Software v1.2 application, “Help” tab link to manual

Related TGL documents:

- YYYY_MM_DD_SAMPLE SUBMISSION FORM_PI_Lastname_Firstname_TGL.xls
- SOP_Qubit_Assay_for_Nucleic_Acid_Quantification.doc
- 16_05_11_Kapa_Hyperprep_lot_tracking_sheet.xls
- 2016-04-22_095434_Dax_plate_cycling_template.eds
- SOP_RT-qPCR_Analysis_in_Excel_Template.xlsx
- SOP_TapeStation_4200_V1.0

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 1 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

2.0 Materials

<i>Workspace</i>		
Post-PCR area		
<i>Equipment and associated consumables</i>		
ThermoFisher	ABI Applied Biosystems ViiA7	
	QuantStudio Real-Time PCR Software v1.2	
	MicroAmp® Optical 384-Well Reaction Plate with Barcode	4309849
	MicroAmp® Fast Optical 96-Well Reaction Plate with Barcode, 0.1 mL	4346906
	MicroAmp® Optical Adhesive Film	4311971
	MicroAmp Adhesive Film Applicator	4333183
BioRad	Sealing Roller	MSR0001
Eppendorf	Eppendorf Centrifuge 5810 R, refrigerated, (use rotor A-4-81)	022625501
	Eppendorf MTP/FLEX Bucket, n=2	022638866
	Eppendorf Centrifuge 5424, non-refrigerated, without rotor, rotary knobs, 230 V/50 – 60 Hz	022620401
Illumina	Illumina Resuspension Buffer (RSP) PN15026770 (-20°C, or 4°C for working stocks)	Component of PN20020599
Roche (KAPA)	Kapa Library Quantification Kit-Illumina platforms ROX Low qPCR Master Mix (includes 6 standards, Primer Mix, qPCR Master Mix), 500 reactions (-20°C)	7960336001 (KK4873)
	KAPA DNA Standards 1-6 (80ul each, 0.0002pM- 20pM), PN7960093001 (-20°C) **may be ordered separately	
	KAPA Primer Mix, 1mL, PN7960093001 (-20°C)	
	KAPA SYBR FAST qPCR Master Mix, 500 reactions (-20°C) (does not contain required primers!)	
Medstore	Nuclease-free water (RT)	W4502-1L
	Tween-20, Biotech grade, Nuclease free, 500ML (RT)	TWN510.500

3.0 General Guidance

- Do not mix and match reagents from multiple kits! Each standard kit is sufficient for 6 RT-qPCR plates. Do not blend standards from different lots!
- When opening a new kit, record all kit lot numbers in TGL master lot tracking sheet. (R:\Lot_tracking_forms\16_05_11_Kapa_Hyperprep_lot_tracking_sheet.docx, associate RT-qPCR excel file with lot letter key (“A”, “B”) for KAPA kit and KAPA standards.
- Before beginning work every day, wipe down all pipets and bench surfaces with peroxide wipes, then 70% ethanol (made from bulk ethanol, 4L) to remove residual detergents.
- Standards and qPCR mix may be briefly vortexed to ensure thorough mixing (exception to rule).
- Thawed standards and aliquoted master mix may be put in the fridge for short term storage or thawing prior to use.

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 2 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

- When loading qPCR plate, all library dilutions and master mix may be kept at room temperature, as chilled reagents will affect pipetting volume.
- PCR master mix should be kept in the dark when not in use to prevent SYBR degradation.
- Liquid in PCR plate may be electrostatic and “jump” to the adhesive covers. Handle carefully.
- Tween-20 is extremely viscous! Aliquot 500ulX10 tubes from Tween-20 PNTWN510.500 stock for general use. Tween-20 H₂O mixes are prone to growing bacteria/fungus; do not store for more than 3 weeks. Always examine solution to confirm no observable growth is present.
- DNA standards are a 10 fold dilution series of a linear 452bp template.
- qPCR primers present in PN7960093001 have the following sequence, primer 1: 5'-AAT GAT ACG GCG ACC ACC GA-3', primer 2: 5'-CAA GCA GAA GAC GGC ATA CGA-3'
- Illumina Resuspension Buffer (RSB) from any Illumina Library Kit may be used for dilution of stock libraries. Aliquot RSP into 1 mL tubes and store in -20°C, working aliquots may be stored at 4°C.

4.0 Procedure

POST-PCR AREA

Preparation for Plate Loading:

1. When opening a new kit, add exactly 1000ul of 10X Primer Premix PN7960093001 to 5mls of provided 2X KAPA SYBR FAST qPCR Master Mix, discard excess primer. Vortex and aliquot 1mL of qPCR-primer mix into six separate 1.5ml microtubes, labeling “KAPA qPCR MM” with date and lot for storage in -20°C and update lot tracking sheet. After each use, indicate freeze thaw with a “dot” on the lid of the tube to track freeze thaw cycles (avoid more than 5 freeze-thaws). Do not blend lots of reagents from different kits!
2. Make a 50ml mix of molecular grade H₂O and 0.05% Tween-20 (25ul of tween-20 in 50mls H₂O). Solution will be used for RT-qPCR 1 in 1000 dilutions. 0.05% Tween-20 solution is stable for 3 weeks at RT.
3. In a dark drawer, thaw 1ml aliquot of premixed RT-qPCR mix (containing primer and master mix, from step 1), standards, and libraries for quantification. Thawing may be done at room temperature if ready, or tubes may be placed in fridge until ready to begin. Briefly vortex “KAPA qPCR MM” and standards, spin down.

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 3 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

Library Dilutions and Standards:

Dilute each stock sequencing library to a library sequencing dilution (MISO DIL) with a concentration 0.5 ng/ul, 50 ul total volume in Illumina Resuspension Buffer (RSP). If stock sequencing library concentration is less than 0.5ng/ul do not dilute; perform RT-qPCR on top stock.

Keep library dilutions as they will be used for sequencing. The “1 in 999ul” dilutions may be discarded after analysis of RT results. See appendix A, reference diagram.

- 1) Each sample library requires 3 separate dilutions, switch tips between dilutions to prevent contamination to top stock. Briefly vortex each stock library and spin down. Dilute each top stock library into 3 replicate dilutions of 1ul library (sequencing dilution) in 999ul of H₂O +0.05% Tween-20 (25 ul of Tween-20 in 50 mls of H₂O). Vortex briefly. Each dilution will be loaded into 2 separate wells on the PCR plate. Therefore each library will use 6 wells of master mix.
- 2) Standards require 18 wells (3 wells per standard, 6 standards).

Plate Loading:


- 1) Each well requires 6 ul of master mix. Briefly vortex master mix and spin down.
- 2) Add 6 ul of master mix to all required wells of PCR plate. You may touch the side of the well with the pipette tip. When all wells are loaded, tap the plate gently to collect pipetted master mix at the bottom of the wells. Seal the plate with MicroAmp adhesive film and plate sealer roller or applicator, and spin at 1000RPM for 1 minute.
- 3) Remove adhesive film without disturbing solution at bottom of wells. If you do, spin the plate down again, and verify no mix was lost.
- 4) Aliquot 4 ul of diluted library (1 ul in 999ul of H₂O + 0.05X Tween-20) or 4ul of standard into PCR well. You may touch the side of the well with the pipette tip. Each standard uses 3 wells, each library dilution requires 2 wells. Change pipette tips between individual library dilutions, and between standards. Seal plate with MicroAmp Adhesive film with plate sealer roller/wedge. Run around the plate margins with the end of the plastic wedge to seal the cover tightly (otherwise evaporation can occur).
- 5) Spin the plate at 1000RPM for 1 minute.
- 6) Wipe the optical surface of the plate with a Kimwipe. Plates may be kept in the fridge for approx. 4 hours before loading, but spin again before loading.

Loading the RT-qPCR Plate in the ViiA7:

- 1) Turn on the Applied Biosystems ABI ViiA7 RT-qPCR machine.
- 2) Ensure that the correct thermal block “Fast 96-Well (0.1mL)” or 384 well block, and PCR plate rack is installed. Blocks are stored in the drawers below the instrument.
- 3) Open the QuantStudio Real-time PCR software which must connect to the instrument.
- 4) Load the plate in the ViiA7 RT-qPCR machine.


SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 4 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

- 5) Open TGL's template file (2016-04-22 095434_Dax_plate_cycling_template.edc) which contains pre-set run conditions for cycling; verify settings below are loaded.

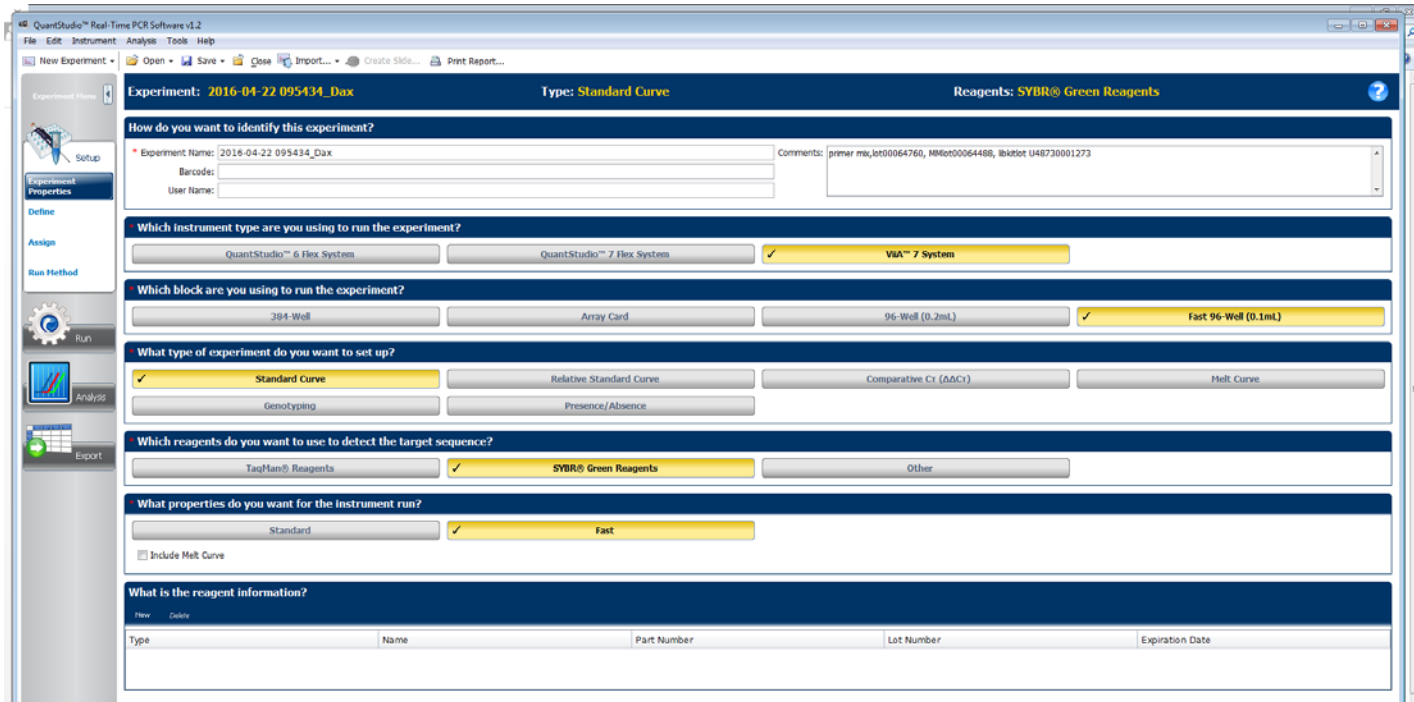

6. Under the run method tab, verify the following settings (see image below)
 - a. Reaction volume =10 ul
 - b. Hold stage= 95C for 5 minutes, data collection off
 - c. PCR Stage= 35 cycles, 95C-30 seconds, 60C- 45 seconds, data collection on

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 5 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

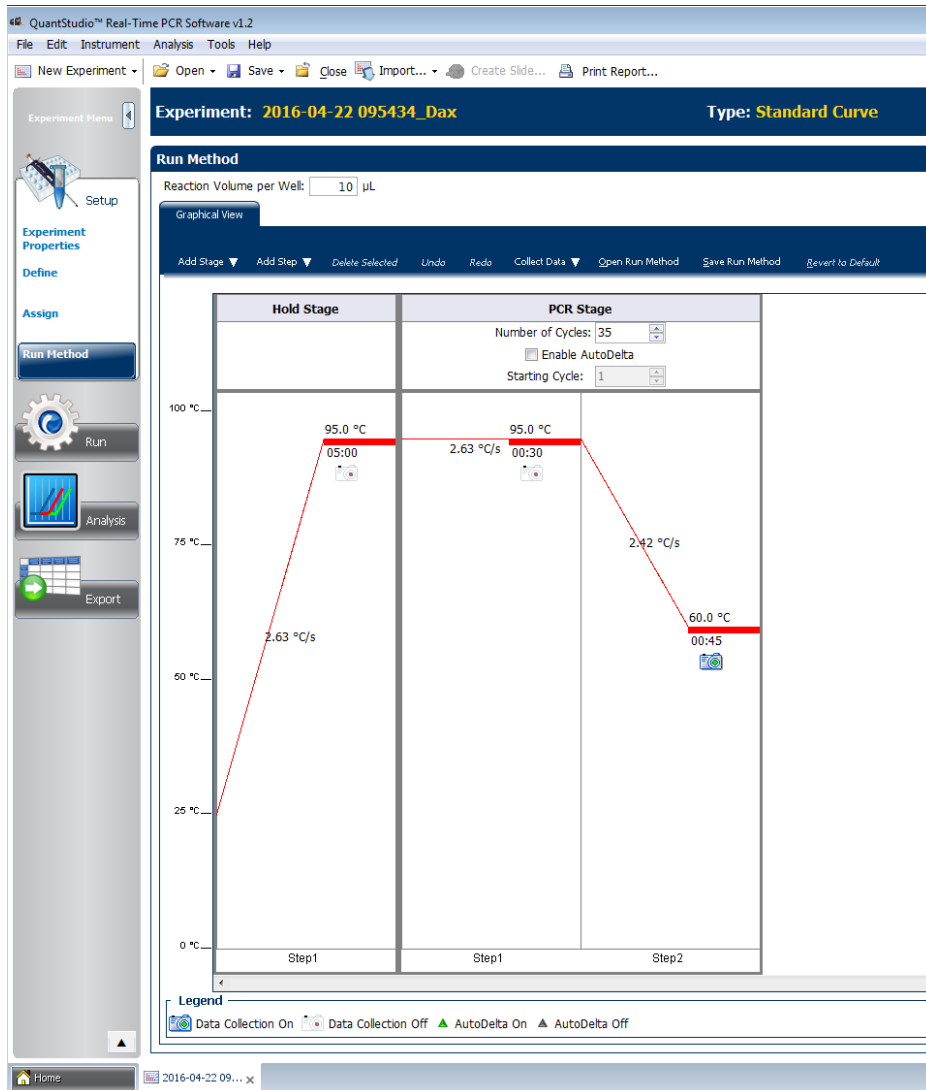


7. Targets may be defined under the “define tab” on the left margin of the window. Individual samples may be indicated here and will export when the run is complete. Ensure all used wells and indicated “targets” use “SYBR” reporter, if SYBR is not indicated sample will not be detected.
8. Assign samples and standards to PCR plate wells. The instrument will not detect/measure a well if it is not indicated with template or standard. Save a new file on the desktop in TGL folder.
9. Ensure instrument is connected (Run status), and open PCR door by pressing open on ViiA7 touch screen to extend loading rack. Place spun, Kim-wiped PCR plate on rack; align plate to the A1 rack placeholder, and close door.
10. Start run (approximately 1.5 hours).

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 6 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

Post Run Analysis and Data Export to Excel:

- Click on “Analysis” menu at top of screen and a new window will open. Click on “Standard curve settings” and verify slope of the line (approx. -3.470, and R2 value approx. 0.999, and efficiency approx. 94%. If CT range of KAPA DNA standards vary significantly, record results in RT analysis excel sheet and flag for review, and or possible repeat of RT-qPCR assay.

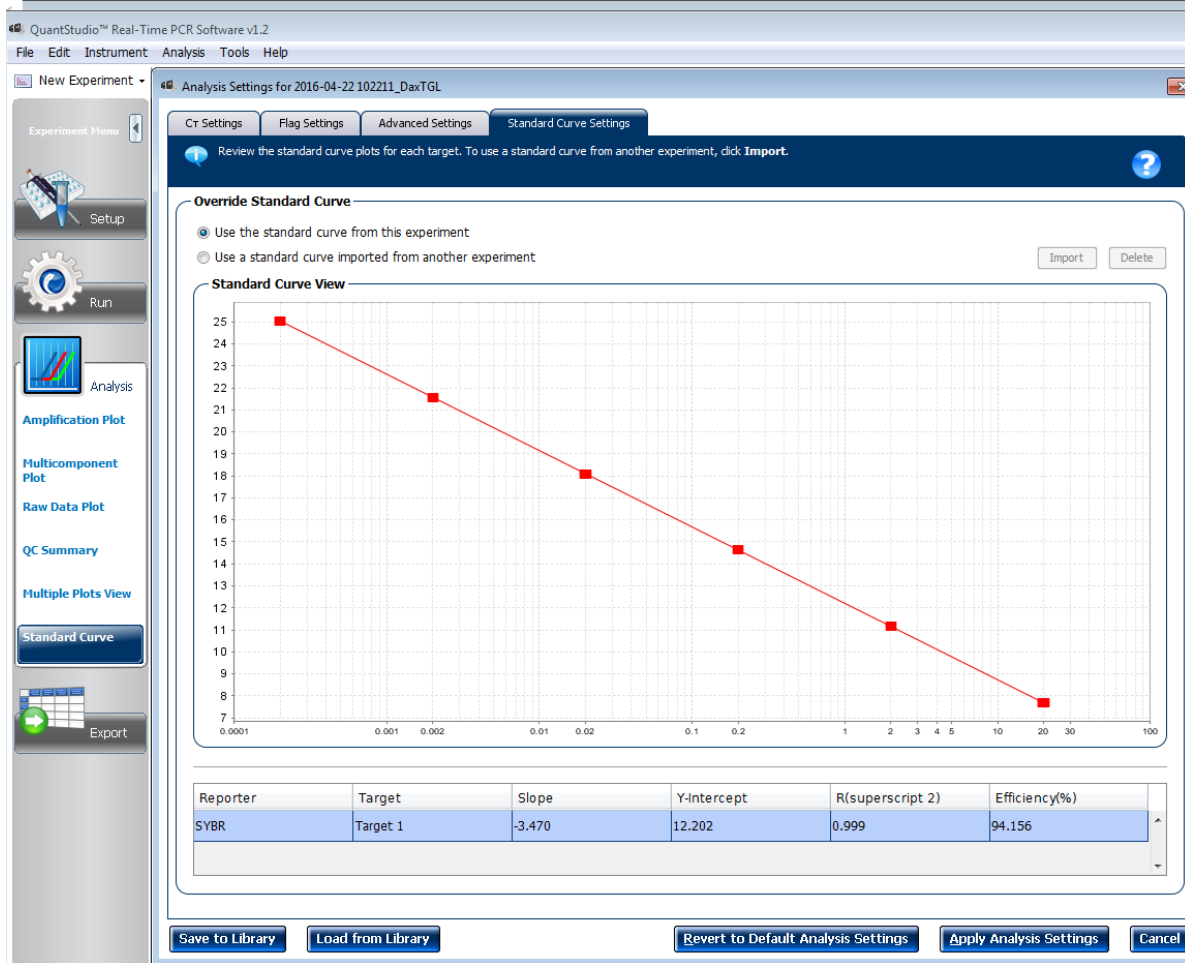

- Email your results (ex: 2016-04-22 102211_DaxTGL.eds) from the instrument PC and store on TGL’s R drive R:\TGL\DNA_QC\RT-qPCR. Do not use USB keys!
- At your desk, open the run “file_name.eds”. Click on the “Export” tab in the left margin. Verify export file location) and export “Results”.
- Use the “SOP_RT-qPCR_Analysis_in_Excel_Template.xlsx” as an example on how to format your excel sheet.
 - Keep a tab “0” with the original export data

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 7 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

- In the run information tab, under notes, record the master mix, standards
- In the calculations tab, take the average of 6 sample wells. CT values should be no more than 0.5 CT apart. Up to two data points may be removed from the average if they fall outside of 0.5 CT from the average.
- Summarize the analysis in the “Summary tab”. Include the TapeStation file name, RT-qPCR file name, and tech signatures. You will need results from the TapeStation to size correct your library using the equation:

$$\text{Final size corrected library concentration (nM)} = \text{pM} * (452 / (\text{average library size from 100-1000 region}))$$

- Record bad results in the summary and note if variability was found! This will be useful if an issue arises with the instrument or reagents that may not currently be apparent.

10. Transfer the summary results to respective sample sheets.

5.0 Quality Control

If CT range of KAPA DNA standards vary significantly (0.5CT average shift for any standard) record results in RT analysis excel sheet and flag for review, and/or possible repeat of RT-qPCR assay. DNA standards should have the following CT +/- 0.5CT:



Table 1: Anticipated CT ranges of KAPA DNA Standards

	Quantity nM	CT
Standard 1	20	7.73
Standard 2	2	10.98
Standard 3	0.20	14.77
Standard 4	0.02	18.02
Standard 5	0.002	21.45
Standard 6	0.0002	24.89

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 8 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

6.0 Appendix A: RT-qPCR Dilution Diagram

Component	Vol (μl)
2x qPCR MasterMix	5
10x Primers	1
1:1,000 subdilution or Standard	4
Total reaction-->	10

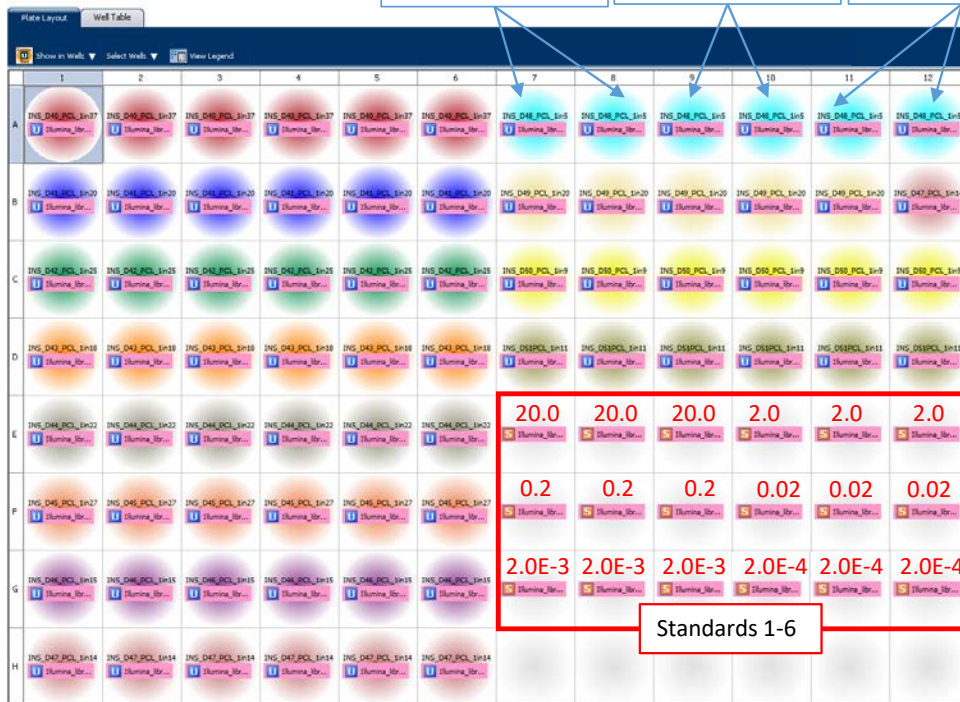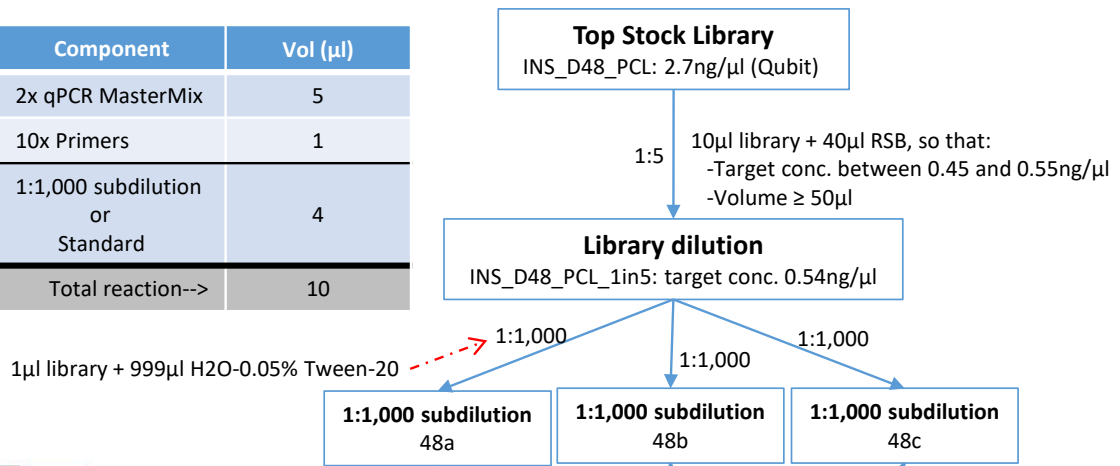

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 9 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	

7.0 Revision History

Version Number	Date (yyyy-mm-dd)	History of change
1.0	2017-05-24	First draft by Dax Torti
1.0.1	2017-05-29	Edited by Alberto Leon
1.0.2	2017-05-29	Kayla Marsh reviewed, in production
1.0.3	2019-07-10	DT- added TapeStation SOP reference. Updated hyperlinks. Added DNA standards reference CT values to Quality Control section.

SOP Title:	KAPA Library Quantification (Illumina Platforms)	
Version:	V1.0.3	Page 10 of 10
Technician approvals		
Approved by:	Dax Torti 190710	
Date:	190710	